


Operation Manual

pH / mV Conductivity / TDS / Salinity Dissolved Oxygen Meter

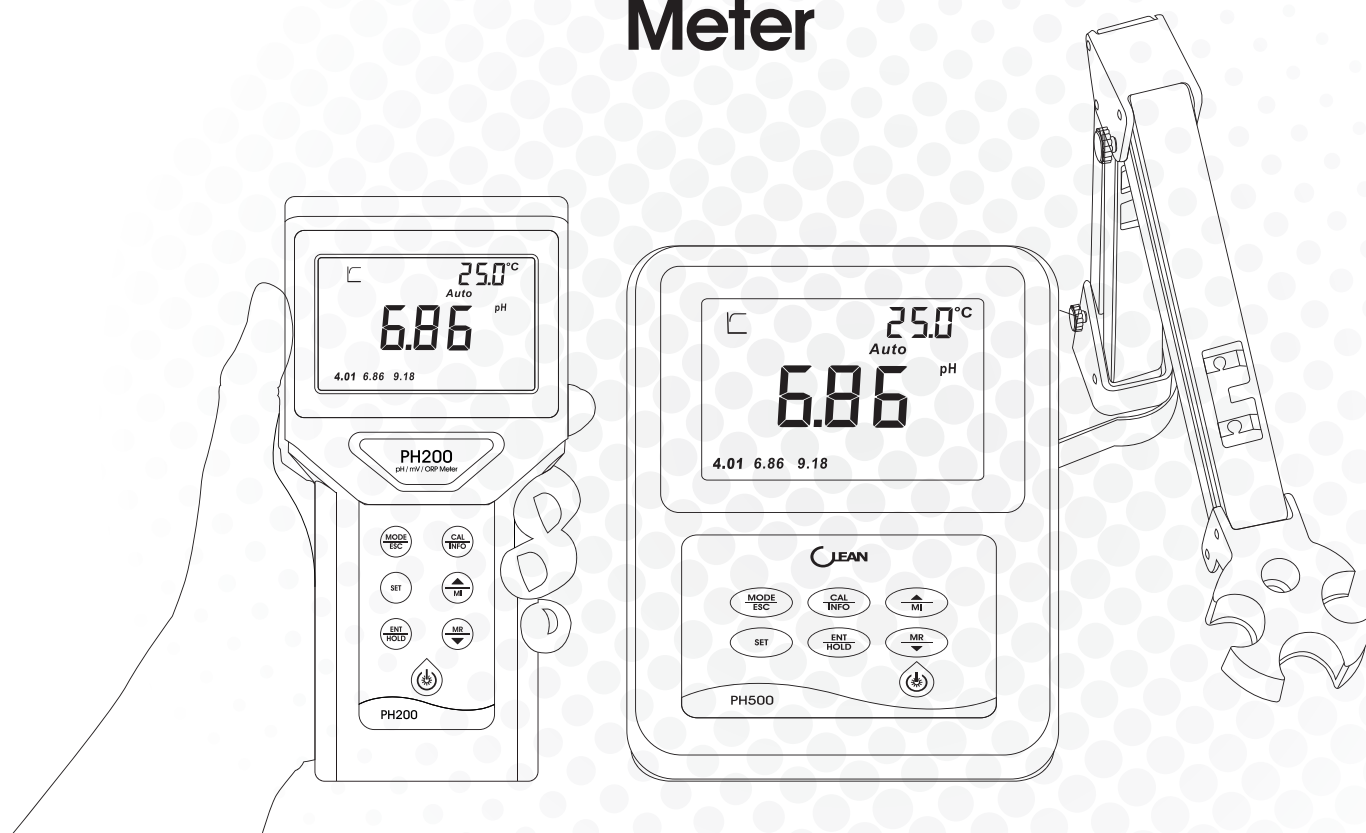

Portable Meter

PH200 / CON200 / DO200

Benchtop Meter

PH500 / CON500 / DO500

PH200 / PH500

CON200 / CON500

DO200 / DO500

TABLE OF CONTENTS

1 Preface	01
2 Basic Operation	02
2.1 Battery	02
2.2 Electrode Installation	02
2.3 Shell	02
2.4 Power on/off and Backlight	02
3 Key Panel	03
3.1 Key instruction	03
3.2 Key function	03
4 Appearance	04
4.1 Display	04
4.2 Display Character Table	05
pH / mV	06
5.1 Technical Specifications	07
5.2 Browse the parameter	07
5.3 pH Set Up	08
5.3.1 P01 pH Standard Solution	08
5.3.2 P02 Manual Temperature setting	09
5.3.3 P03 Temperature Offset setting	10
5.3.4 P04 Auto Lock Function setting	11
5.3.5 P05 Auto Power Off Setting	12
5.3.6 P06 Clearing Memory Function	12
5.3.7 P07 Reverting to Factory Default Setting	13
5.4 Calibration Mode	14
5.5 Measurement Mode	15
5.6 Storage of Data	16
5.7 Calibration Parameter and Setting Parameter	17
5.8 Appendix	18
Conductivity / TDS / Salinity	19
6.1 Technical Specifications	20
6.2 Browse the parameter	21
6.3 Set Up	22
6.3.1 P01 Cell Constant (k)	22
6.3.2 P02 Temperature Compensation Base	23
6.3.3 P03 Conductivity Temperature Factor	23
6.3.4 P04 TDS Factor	24

6.3.5 P05 Temperature modified Factor	24
6.3.6 P06 Auto Lock Function Setting	25
6.3.7 P07 Auto Power Off Setting	25
6.3.8 P08 Clearing Memory Function	26
6.3.9 P09 Reverting to Factory Default Setting	26
6.4 Setting Parameter	27
6.5 Calibration	28
6.6 Appendix	29
6.6.1 Conductivity sensor measuring range	29
6.6.2 Factory Defaults	29
Dissolved Oxygen	30
7.1 Technical Specifications	31
7.2 Browse the parameter	32
7.3 Set Up	32
7.3.1 P01 Atmospheric Pressure	32
7.3.2 P02 Salinity Compensation	33
7.3.3 P03 Temperature Offset	33
7.3.4 P04 Auto Lock Function Setting	34
7.3.5 P05 Auto Power Off Setting	34
7.3.6 P06 Clearing Memory Function	35
7.3.7 P07 Reverting to Factory Default Setting	36
7.4 Calibration	37
7.5 Measurement Mode	38
7.6 Storage of Data	39
7.6.1 Storage of measuring data	39
7.6.2 Browse on data storage	39
7.6.3 Setting parameter	40
7.7 APPENDIX	41
7.7.1 The corresponding Average Atmospheric Pressure with the Altitude	41
7.7.2 Water saturated dissolved oxygen in different temperature	42
7.7.3 Factory defaults	42
8 GENERAL INFORMATION	43
8.1 Warranty	43
8.2 Return Of Malfunction Instruments	43
8.3 Guidelines Or Returning Unit For Repair	43

1 Preface

Thank you for selecting CLEAN meter.

Although the meter use advanced technology and meet the requirements of current safety rules, improper use can still threaten the safety of users, and / or cause harmful influences to factory and other equipments. Therefore, before using the meter, relevant person must read and understand contents of this operation manual.

Besides the reasonable prices, Clean meters also have the following advantages:

- Simple operation. This operation manual provides you clear and easy operation guidance.
- Humanized design, Comfortable holding.
- Multiple accessories available, such as sensors, buffer solution etc.

	Portable Meter	Benchtop meter
PH	<input type="checkbox"/> PH200	<input type="checkbox"/> PH500
Conductivity / TDS / Salinity	<input type="checkbox"/> CON200	<input type="checkbox"/> CON500
Dissolved Oxygen	<input type="checkbox"/> DO200	<input type="checkbox"/> DO500

Safety



- Never expose the meter in the explosive environment! Part of the shell case is not airtight and the invasion of spark or the corrosion caused by the invasive gas may cause explosion.



- Please follow the operation manual and the laboratory safety regulations when use the chemicals and solutions.

Safety precautions



- Never split the shell case.
- Only the OEM maintenance staff is allowed to maintain the meter.
- Following environment effect should be avoided:
 - Violent vibration
 - Expose to sunlight for a long time
 - Atmospheric humidity more than 95%
 - Corrosive gas
 - Ambient temperature below than -10°C or more than 60°C
 - Strong electric or magnetic field

2 Basic Operation


2.1 Battery — (200 Series)

Use a coin to remove the battery cover by following the “-” direction.

Insert 2*1.5V 7AAA batteries in the battery compartment with the correct Positive and Negative poles.

Tight the battery cover by following the “+” direction.


PH200 / CON200 / DO200

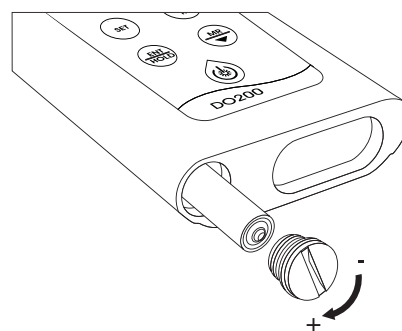


2.2 Electrode Installation

You can refer to the Electrode operation manual for use and maintenance. The direction of the red dot marker on the Electrode and the salient point on the unit should be the same when install the electrode to the unit. Remove the electrode by pulling out the clamp on the electrode.

PH200 / PH500

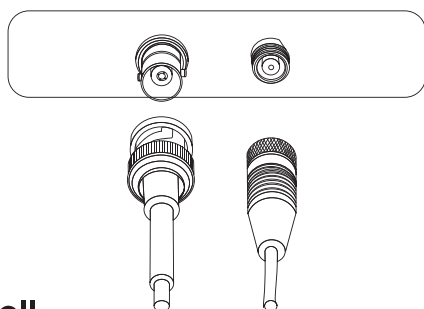

CON200 / DO200 / CON500 / DO500

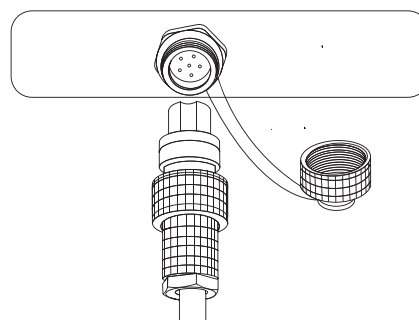

2.3 Shell

This series of meter is portable . If the meter accidentally touches water or other fluids, please wipe it clean immediately. The meter function is not affected usually. But if affected and cannot use, please contact our company or CLEAN local customer service center. They will be very glad to help you.

2.4 Power on/off and Backlight

Power on

Short press power on/off key to start the meter. The whole screen displayed within 1 second.

Power off

Long press the power on/off key to power off the meter.

Auto power off in 10 minutes if without any operation.

LCD backlight

Short press power on/off key to control if turn on the backlight or not.

3 Key Panel








3.1 Key instruction

Short press: Short press means release the key once after pressing. (If there is no mark out below, default it as Short press)

Long press: Long press means press the key for 3 seconds and release.


Hold: Hold means not release the key, used in adjusting data and accelerate after a certain time. Not release the key until the value is set as needed.

3.2 Key function:

Key	Description
	Power on ON/OFF backlight Long press to Power off
	Mode switch Exit from current mode operation
	Data setting
	Enter into calibration mode Long press to enter into parameter browsing
	Freeze or unlock the displayed value Press to unlock in measurement mode
	Store the displayed value into memory Increment values or scroll through the next options available Press to browse the saved data in measurement mode
	Recall stored values from the memory Decrement values or scroll through the next options available Press to browse the saved data in measurement mode

4 Appearance

4.1 Display

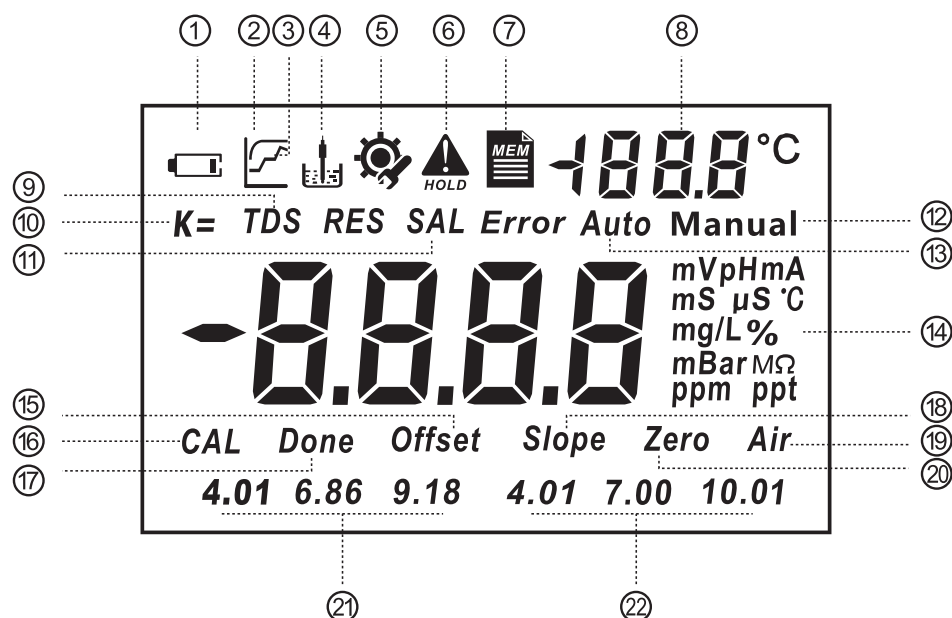

- 1 Battery power
- 1 Measuring Status - Calculating
- 2 Measuring Status - Stable Value
- 3 Electrode inserted display
- 4 Setup display
- 6 Auto lock
- 7 Memory
- 8 188.8°C Temperature
- 9 TDS - Total dissolved solids
- 10 k= - Cell Constant
- 11 SAL - Salinity
- 12 Manual - Manual Temperature Compensation
- 13 Auto - Automatic Temperature Compensation
- 14 mV, pH, mA, °C, %, mS, μS, °C, mg/L, ppm, ppt, g/L, mBar - Unit of Measurement
- 15 Offset - Electrode Offset
- 16 CAL - Calibration
- 17 Done - Calibration Done
- 18 Slope - Electrode Slope
- 19 Air - Air Calibration
- 20 Zero - Zero Calibration
- 21 4.01, 6.86, 9.18 Calibrated Points - NIST Buffer Standard
- 22 4.01, 7.00, 10.01 Calibrated Points - USA Buffer Standard

4.2 Display Character Table

SEN	Sensor	CODE	Password Setting
t.S	Manual Temperature setting	DEF	Factory Defaults
t.C	Automatic Temperature Compensation	SL1	Sensor Slope 1
Auto	Automatic Temperature Compensation	SL2	Sensor Slope 2
Manual	Manual Temperature Compensation	SAVE	Save Data
Slope	Electrode Slope	Err	Error
Hd	Autolock	ON	On
ADF	Auto Off	OFF	Off
NONE	no stored data	NO	No
CLr	Clear data	YES	Yes
d-	place of the stored data	Over	Temperature value Over
Offset	Sensor Offset	Under	Temperature value Under
P-	Menu item	FULL	Full Data Storage
		Over	Measuring Value Over
buf	Buffer Solution	Under	Measuring Value Under
NIST	NIST Standard		
USA	USA Standard	NA	nA output of electrode
		SLP	Sensor slope
t.dS	TDS	PrE	Atmospheric pressure
SAL	Salinity	SAL	Salinity
t.rF	Temperature Base	OSP	%
COE	Coefficient	OC	mg/L
OFFS	Offset	OPr	DO electrode voltage

pH / mV

PH200 Portable Meter

PH500 Benchtop Meter

2014 New Model! Brief and exquisite design, space saving, easy calibration with calibrated points displayed, optimum accuracy, simple operation comes with high luminant backlight. It is your reliable partner for routine applications in laboratories, productions plants and schools.

PH200 / PH500

- All-weather precise, Comfortable holding, Easy carrying , Simple Operation
- 65*40mm, large LCD with backlight for easy meter information reading
- IP67 rated, dustproof and waterproof, Floats on water - **[PH200]**
- Up to 3 points auto-buffer-recognition: Zero offset, Slope of Acid/Alkali segment, ensure accurate measuring results for the full range
- One key to check through all the settings, including: Zero offset, Slope of Acid/Alkali segment, and all the settings
- Auto lock function
- 50 data sets memory stores and recalls
- Auto Power off saves battery after 10-mins non-use
- 2*1.5V 7AAA battery, long battery life - **[PH200]**
- CP337 Operating Carrying Pouch - **[PH200]**
- Economical, easy to operate, cost saving
- Detachable Electrode Stand organizes multiple electrodes neatly, easy installing at either left or right side and holds them firmly in place - **[PH500]**

5.1 Technical Specifications

		PH200	PH500
pH	Range	-2.00 ~ 16.00 pH	
	Resolution	0.01 pH	
	Accuracy	±0.2% F.S.	
ORP	Range	-2000~2000 mV	
	Resolution	1 mV	
	Accuracy	± 0.2% F.S.	
Temperature	Range	-10.0 °C~110.0 °C	
	Resolution	0.1 °C	
	Accuracy	± 0.2 °C	
Power	Power supply	2*1.5V 7AAA batteries	Transformer 100 ~ 240V
	Working time	>500 hrs	—
pH Buffer Kinds	NIST	4.00、6.86、9.18	
	USA	4.01、7.00、10.01	
Others	Environment	-5 °C~60 °C, Relative humidity<90%	
	Memory	50 sets	
Dimensions		94*190*35mm (W*L*H)	140*210*35mm (W*L*H)
Weight		250g	650g

5.2 Browse the parameter


Long press  for 3 seconds in measurement mode.

Then press  or  to browse the parameters.

1. Zero offset
2. Slope
3. High/Low segment of the slope
4. Buffer kinds
5. Temperature compensation
6. Hold, Auto Lock
7. Auto power off
8. Battery power - **[PH200]**

5.3 pH Set Up

5.3.1 P01 pH Standard Solution

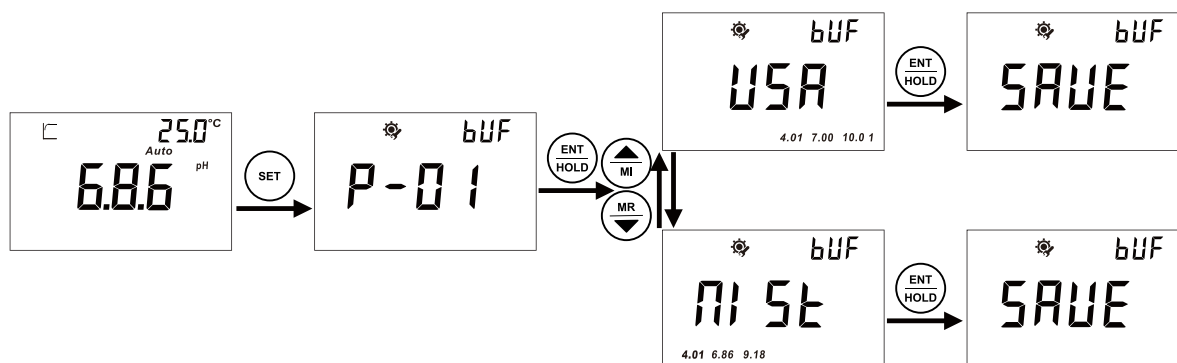

- After power-up, the unit will enter the measurement mode. Factory default is pH measuring mode. You can switch between the pH measurement mode and the ORP measurement mode by pressing the Key.
- Press to enter SET up step; Press to enter the Electrode Calibration mode.
- In P-01, you can select a group of buffer solutions as a standard: USA (4.01、7.00、10.01) or NIST (4.01、6.86、9.18).
- Please refer to above description steps to set up P-01.
- You can go to next parameter setting by pressing , or pressing key to quit and go back to measurement mode.
Factory default: USA

Note:

1. During setting, you can press key to quit setting whenever necessary.
2. During setting, you can adjust anything twinkling by pressing . If it is the data twinkling, you can speedy adjust the data by pressing .

5.3.2 P02 Manual Temperature setting










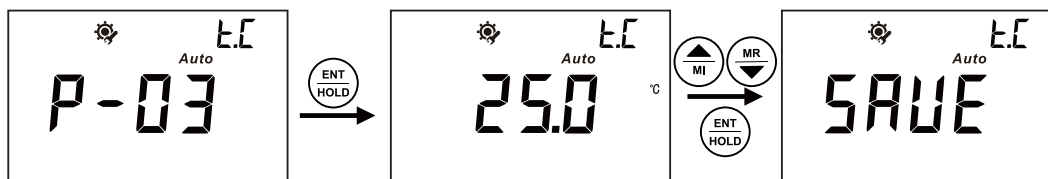

- After entering P-02, you can set the manual temperature by pressing . Long press for speedy setting. Confirm your setting by pressing .
- Manual Temperature setting range: 0°C-100°C.
- Please refer to above description steps to set up P-02.
- You can go to next parameter setting by pressing , or pressing key to quit and go back to measurement mode.

Note:

1. Please make sure the Temperature Electrode is UNPLUGGED when manual Temperature compensation. Otherwise, the unit will display “Err” and go back to the set up menu for fresh operation. The “Manual” is displayed beneath the main screen normally.
2. It is not working to unplug the Temperature Electrode while setting. You have to set again after removing the electrode and quit the current mode.

5.3.3 P03 Temperature Offset setting











- After entering P-03, you can set the temperature to the appropriate value by pressing   according to the actual temperature. Long press   for speedy setting. Confirm your setting by pressing .
- Temperature Offset range: +10°C.
- Please refer to above description steps to set up P-03.
- You can go to next parameter setting by pressing  , or pressing  key to quit and go back to measurement mode.

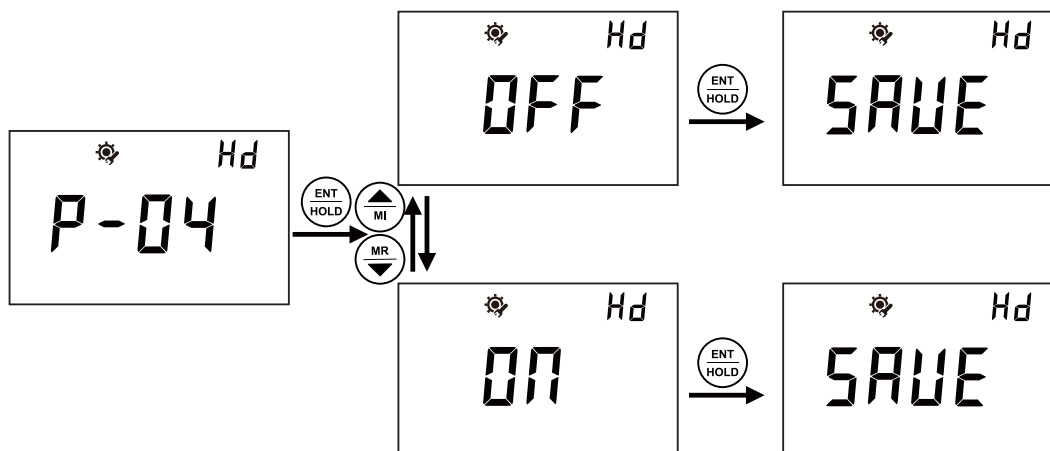Note:


1. Please make sure the Temperature Electrode is PLUGGED when Temperature Offset setting. Otherwise, the unit will display "Err" and go back to the set up menu for fresh operation. The "Auto" will be displayed beneath the main screen the screen normally.
2. It is not working to plug the Temperature Electrode while setting. You have to set again after plugging in the electrode and quit the current mode.

5.3.4 P04 Auto Lock Function setting




- After entering P-04, you can select ON or OFF the Auto Lock Function by pressing  . If the Auto Lock Function is ON, the meter can lock the measured value after the reading had stabilized and display . Press  to unlock.
- Please refer to above description steps to set up P-04.
- You can go to next parameter setting by pressing  , or pressing  key to quit and go back to measurement mode.

5.3.5 P05 Auto Power Off Setting

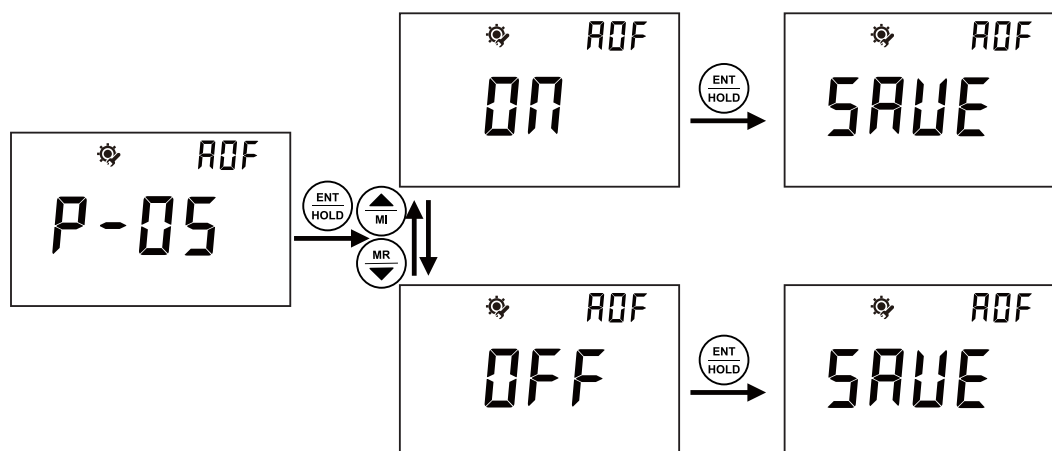


- After entering P-05, you can select ON or OFF the Auto Power Off by pressing . If the Auto Power Off is ON, the meter will power off if no operation in 10mins.
- Please refer to above description steps to set up P-05.
- You can go to next parameter setting by pressing , or pressing key to quit and go back to measurement mode.

5.3.6 P06 Clearing Memory Function

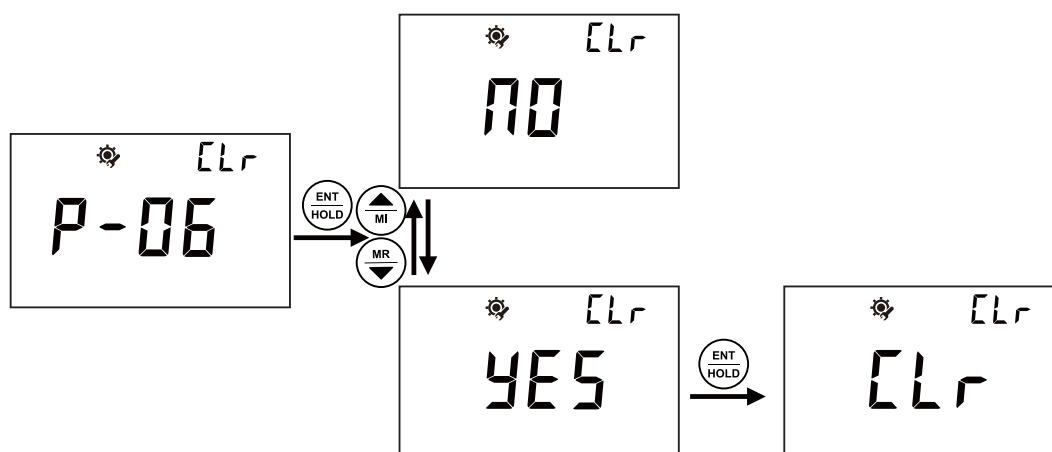

- After entering P-06, you can select NO or YES to not clear or clear the data by pressing . If select YES, the meter will clear all the saved data.
- Please refer to above description steps to set up P-06.
- You can go to next parameter setting by pressing , or pressing key to quit and go back to measurement mode.

5.3.7 P07 Reverting to Factory Default Setting

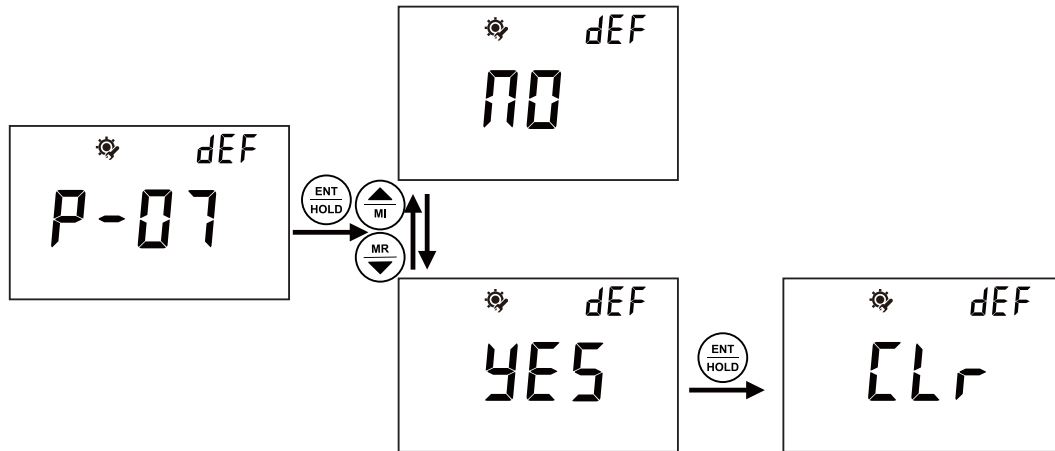


- After entering P-07, you can select NO or YES to not reverting or reverting to factory settings by pressing . If select YES, the meter will revert all the settings to factory defaults and all the settings will lost forever. The unit will restart at the same time.
- Please refer to above description steps to set up P-07.
- You can go to next parameter setting by pressing , or pressing key to quit and go back to measurement mode.

Note:

1. During setting, you can press key to quit setting whenever necessary.
2. During setting, you can adjust anything twinkling by pressing . If it is the data twinkling, you can speedy adjust the data by pressing .

5.4 Calibration Mode

1 pH Calibration

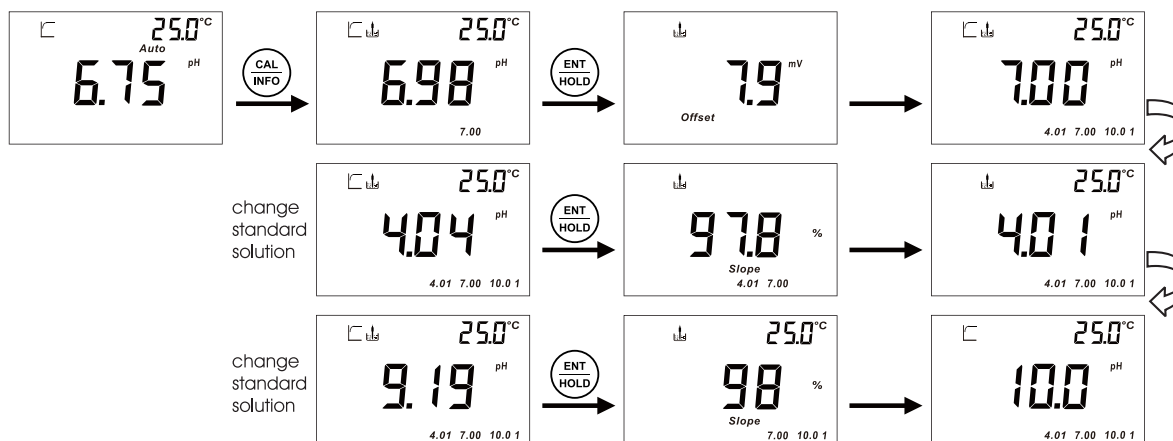

2. Insert the electrode in the pre-set standard solution. You can refer to P-01 explanation on the standard solution setting.

3. Press **CAL INFO** and wait the measuring value to be stabilized. Then press **ENT HOLD** to confirm. The offset slope will be shown after the calibration.


Note: System defaults first point calibration is 6.86pH or 7.00pH.


After the first point calibration, the second point calibration will begin automatically. Then the third point calibration will be continued. The meter will back to measurement mode after all the calibration is done.

Note:




You can press **MODE ESC** back to the measurement mode during calibration and all the calibration point is saved. When the measured signal is out-ranged, the sign "Err" will be showed, possibly the aging electrode or the polluted standard solutions lead to this.

5.5 Measurement Mode

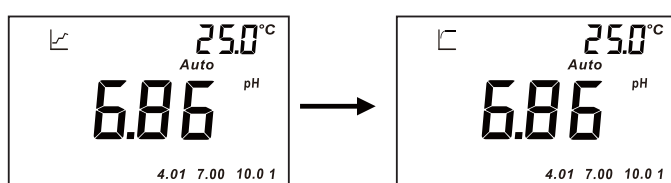1. You can first set the meter per your request. Otherwise, the setting will be factory defaults.
2. Use tap water or distilled water to clean the pH and the temperature electrode, removing the adhesive impurities on the surface of electrode.
3. After power on, press  to select the operation mode (pH or mV).
4. Immerse the pH and the temperature electrode in the solution and stir gently. The measuring mark will twinkle. You can read the value once the data is stabilized and shows the stable mark.






Note:

1. When the screen shows “Auto” means the temperature sensor is working and shows the actual temperature upper right corner. You can refer to the “Temperature Offset setting” if a revision on the temperature is required.
2. If the “Manual” is shown below the screen means the temperature electrode is disconnected and is manual temperature compensation. Factory default is 25°C. You can refer to P-02 “Manual temperature setting” to adjust the temperature.
3. The measuring value is stabilized when the screen showed  , The  is showed if the Auto Lock Function is on. Press  to unlock.

5.6 Storage of Data


5.6.1. Storage of measuring data







1. If you need to save the measuring data once it is stabilized, press  , the Lot no. will be shown on the upper right of the screen. Press  to save the data. You can also press  to not save. The meter will automatically go back to measurement mode after the data saved. The unit at most can save 50 sets data. The screen will show "OVER" if more than 50 sets reminds you the memory card is full.

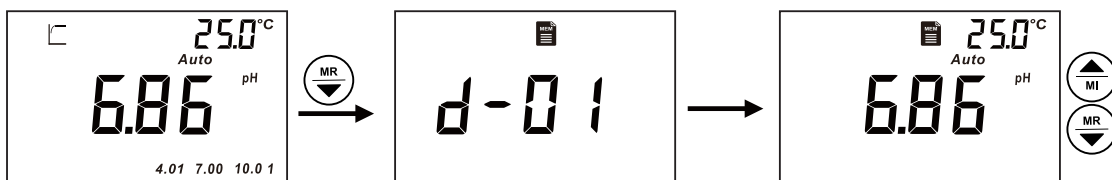2. Enter into setting mode P-06 if you need to delete the saved data forever. You can refer to Clearing memory function.


5.6.2. Browse on data storage



1. Press  to enter into the browse on the data storage , the Lot no. will be shown and data twinkling. Press   to check all the saved data. Press  to go back to measuring mode. If showed "NONE" means no data is saved yet.

2. If you quit and re-entering during checking the data, the meter will freeze in the last set of data.

5.7 Calibration Parameter and Setting Parameter

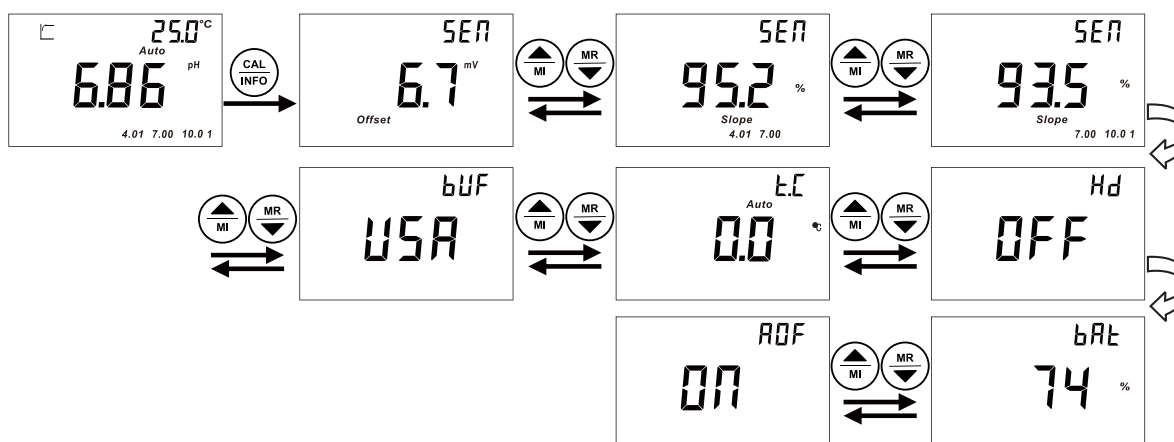

1. Long press to enter into Setting parameter when in measurement mode. Press to check through all the Calibration parameters and Setting parameter.
2. Press to quit and go back to measurement mode.

5.8 Appendix

Factory defaults

Item	Corresponding interface	Factory Defaults
Standard Solution	P-01	USA: 10.01,7.00,4.01
Temperature	P-02	Manual Temperature Compensation 25 °C
Auto Lock	P-04	OFF
Auto power off	P-05	ON
Backlight		OFF

Conductivity / TDS / Salinity

CON200 Portable Meter

CON500 Benchtop Meter

2014 New Model! Compact and humanized design, easy and quick calibration, optimum accuracy in Conductivity, TDS and Salinity measurements, easy operation comes with high luminant backlight make the instrument an ideal research partner in laboratories, productions plants and schools.

- All-weather precise, Comfortable holding, Easy carrying , Simple Operation
- 65*40mm, large LCD with backlight for easy meter information reading
- IP67 rated, dustproof and waterproof, Floats on water - **[CON200]**
- Measuring range: 0.000 us/cm-400.0 ms/cm, Automatic range switching
- Unit display: us/cm;ms/cm,TDS(mg/L), Sal((mg/L),°C
- One key to check through all the settings, including: cell constant, slope and all the settings
- Auto lock function
- 50 data sets memory stores and recalls
- Auto Power off saves battery after 10-mins non-use
- 2*1.5V 7AAA battery, long battery life - **[CON200]**
- CP337 Operating Carrying Pouch - **[CON200]**
- Economical, easy to operate, cost saving
- Detachable Electrode Stand organizes multiple electrodes neatly, easy installing at either left or right side and holds them firmly in place - **[CON500]**

6.1 Technical Specifications

		CON200	CON500
Conductivity	Range	0.000 $\mu\text{S}/\text{cm}$ ~ 400.0 mS/cm	
	Resolution	0.001 $\mu\text{S}/\text{cm}$ ~ 0.1 mS/cm	
	Accuracy	$\pm 0.5\%$ FS	
TDS	Range	0.000 mg/L ~ 400.0 g/L	
	Resolution	0.001 mg/L ~ 0.1 g/L	
	Accuracy	$\pm 0.5\%$ FS	
Salinity		0.0 ~ 260.0 g/L	
		0.1 g/L	
		$\pm 0.5\%$ FS	
		0.65	
Temperature	Temperature	- 10.0 $^{\circ}\text{C}$ ~ 110.0 $^{\circ}\text{C}$	
	Resolution	0.1 $^{\circ}\text{C}$	
	Accuracy	± 0.2 $^{\circ}\text{C}$	
Power	Power supply	2*1.5V 7AAA batteries	Transformer 100 ~ 240V
	Working time	>500 hrs	—
Others	Environment	-5 $^{\circ}\text{C}$ ~ 60 $^{\circ}\text{C}$, Relative humidity<90%	
	Memory	50 sets	
Dimensions		94*190*35mm (W*L*H)	140*210*35mm (W*L*H)
Weight		250g	650g

6.2 Browse the parameter


Long press  for 3 seconds in measurement mode.

Then press  or  to browse the parameters.

1. Cell constant
2. Slope
3. Base temperature
4. Temperature change rate
5. TDS
6. TC Manual Temperature compensation
7. Hold, Auto Lock
8. Auto power off
9. Battery power - **[CON200]**

6.3 Set Up

6.3.1 P01 Cell Constant (k)

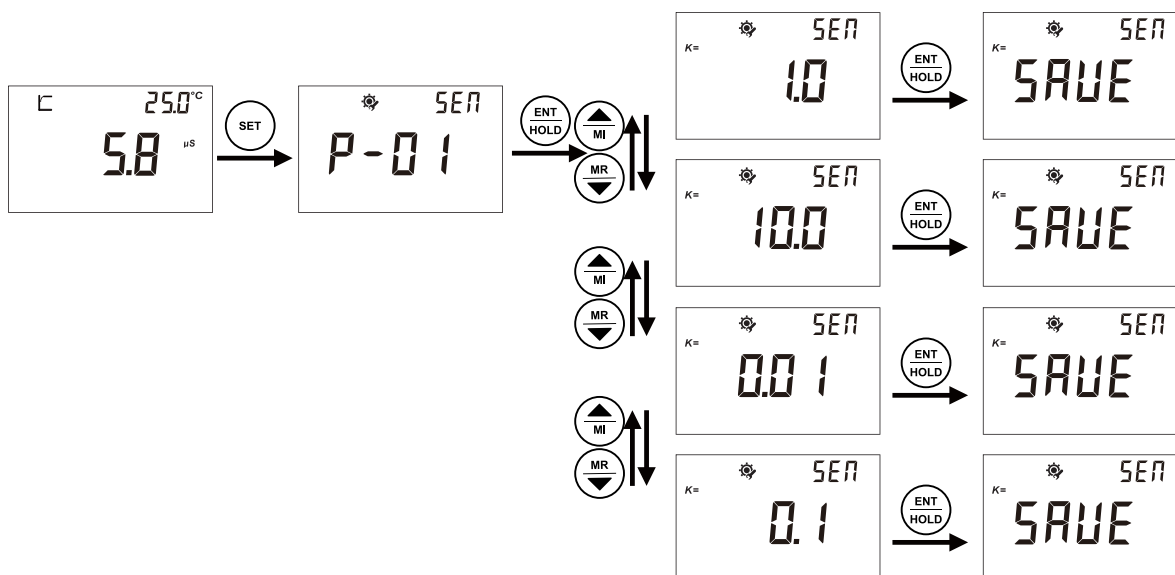

- The unit will enter into measurement mode after power on. Factory default is Conductivity measurement mode. Press **MODE/ESC** to switch between Conductivity measurement mode and Total Dissolved Solids measurement mode. Press **ENT/HOLD** to enter into parameter setting mode. Press **CAL/INFO** to enter into electrode calibration mode.

In P-01, you can set up cell constant k of your sensor for measurement application.

- Please refer to above description steps to set up P-01.
- You can go to next parameter setting by pressing **MI** **MR**, or press **MODE/ESC** key to quit and go back to measurement mode.

Measurement range:

Cell constant k=1.0, For Middle range;

Cell constant k=10, For Wide range (More than 40 mS/cm or 10ppt);

Cell constant k=0.1, For Middle range (Below 40 µS/cm or 10ppm);

Cell constant k=0.01, For Middle range (Below 10 µS/cm) ;




Factory default: Cell constant k=1.0.

Note:

- During setting, you can press **MODE/ESC** key to quit setting whenever necessary.
- During setting, you can adjust anything twinkling by pressing **MI** **MR**. If it is the data twinkling, you can speedy adjust the data by pressing **MI** **MR**.




6.3.2 P02 Temperature Compensation Base



- In P-02, you can set system base value for conductivity measurement.
- Please refer to above description steps to set up P-02.
- Range of Temperature Compensation Base: 15.0-35.0°C.
- You can go to next parameter setting by pressing   , or press  key to quit and go back to measurement mode.




6.3.3 P03 Conductivity Temperature Factor




- In P-03, you can set Temperature coefficient for conductivity.
- Please refer to above description steps to set up P-03.
- Range of Conductivity Temperature Factor: 0.00-10.00%.
- You can go to next parameter setting by pressing   , or press  key to quit and go back to measurement mode.




6.3.4 P04 TDS Factor




- In P-04, you can set TDS factor for your application.
- Please refer to above description steps to set up P-04.
- TDS Factor range: 0.40-1.00.
- You can go to next parameter setting by pressing   , or press  key to quit and go back to measurement mode.

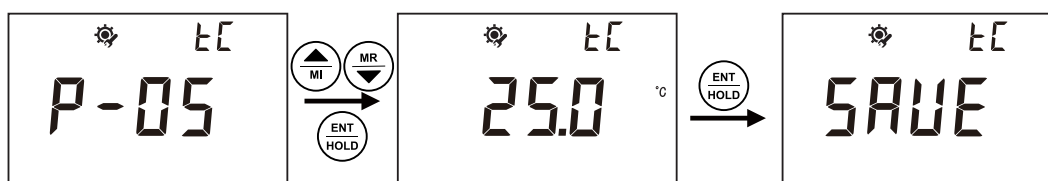6.3.5 P05 Temperature modified Factor




- In P-05, you can set Temperature modified factor for your application.
- Please refer to above description steps to set up P-05.
- Range of Temperature modified Factor: +10°C.
- You can go to next parameter setting by pressing   , or press  key to quit and go back to measurement mode.

6.3.6 P06 Auto Lock Function Setting

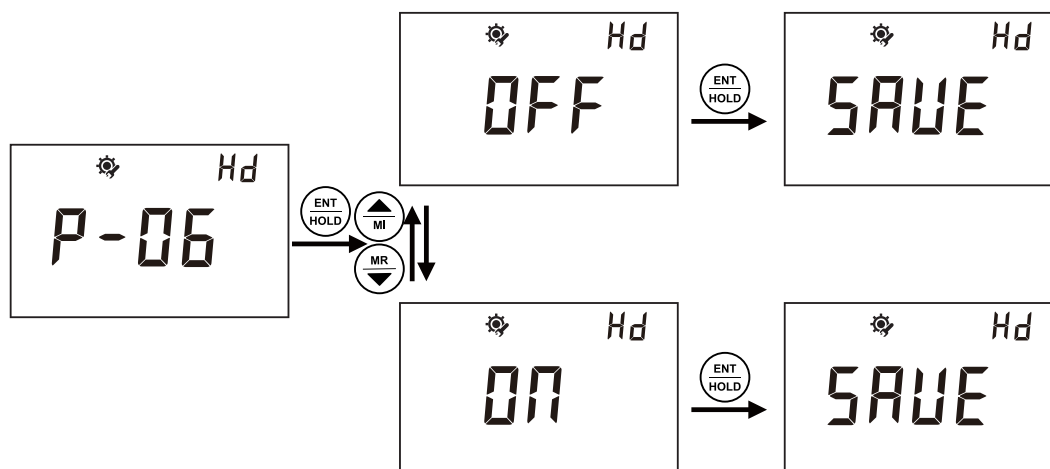


- In P-06, you can set the Auto Lock Function.
- Please refer to above description steps to set up P-06.
- If the Auto Lock Function is ON, the meter can lock the measured value after the reading had stabilized. Press to unlock.
- You can go to next parameter setting by pressing , or press key to quit and go back to measurement mode.

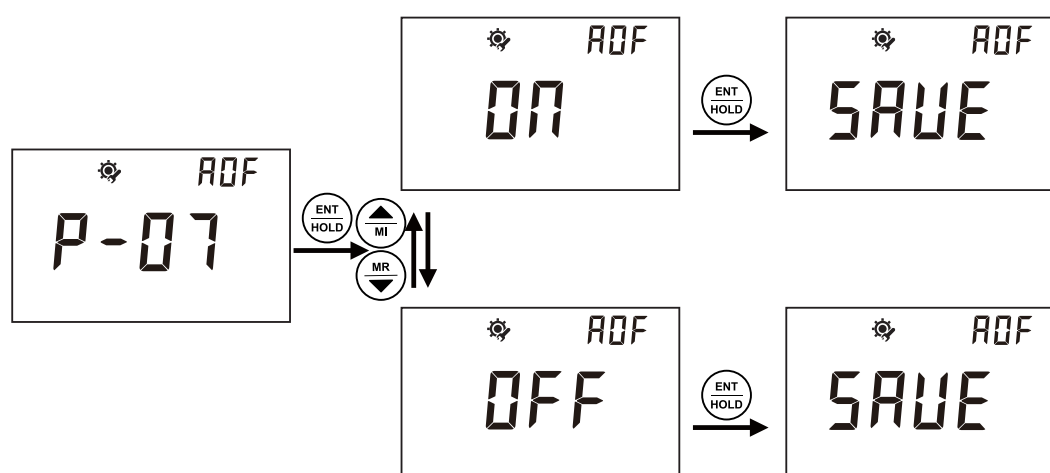


6.3.7 P07 Auto Power Off Setting




- In P-07, you can set the Auto Power Off Function.
- Please refer to above description steps to set up P-07.
- If the Auto Power Off is ON, the meter will power off if no operation in 10mins.
- You can go to next parameter setting by pressing , or press key to quit and go back to measurement mode.

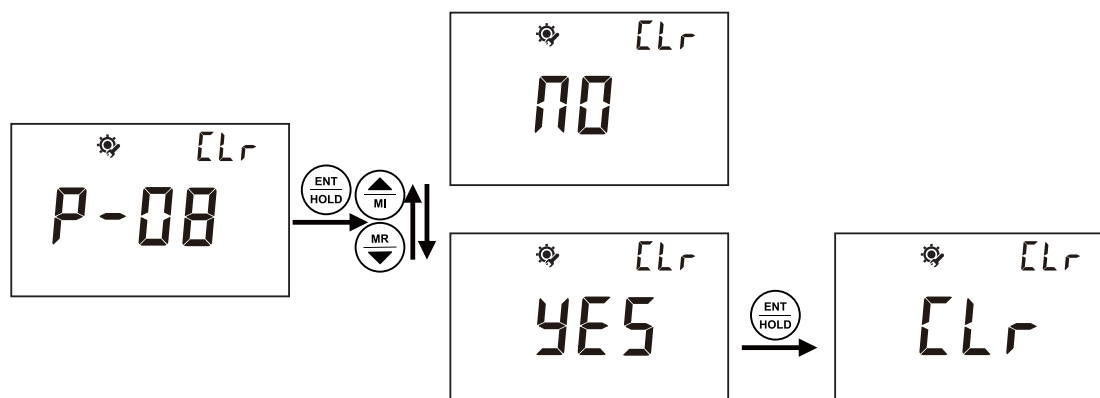




6.3.8 P08 Clearing Memory Function

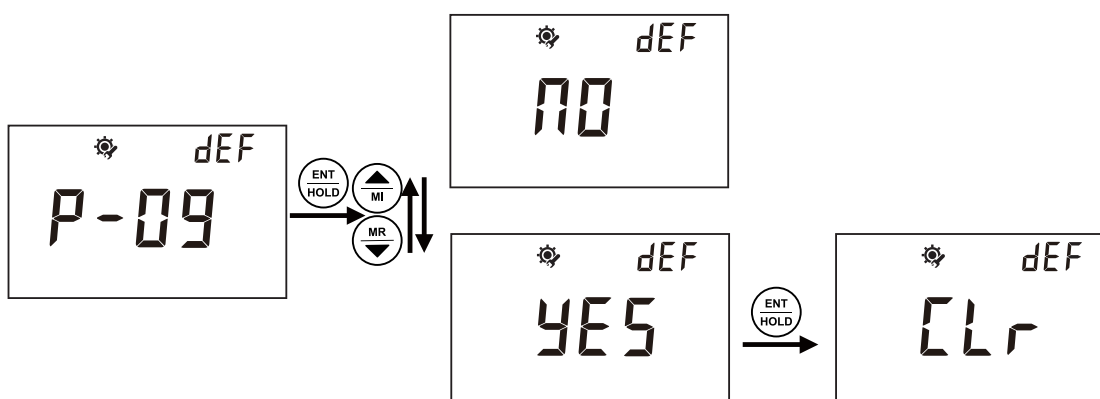


- In P-08, you can set the Clearing Memory Function.
- Please refer to above description steps to set up P-08.
- If select YES, the meter will clear all the saved data.
- You can go to next parameter setting by pressing  , or press  key to quit and go back to measurement mode.





6.3.9 P09 Reverting to Factory Default Setting

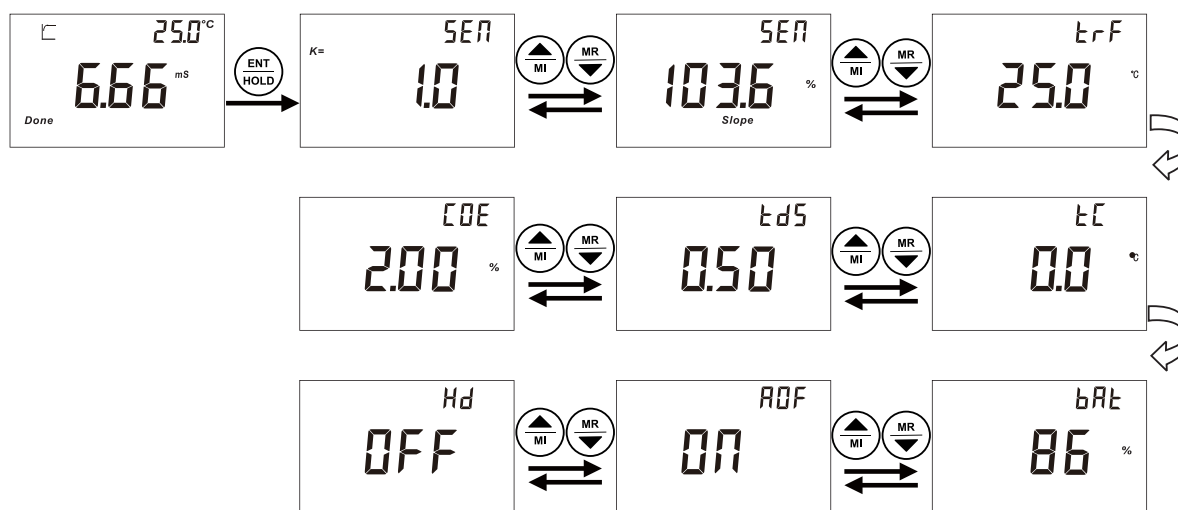

- In P-09, you can select NO or YES to not reverting or reverting to factory settings by pressing  .
- Please refer to above description steps to set up P-09.
- If select YES, the meter will revert all the settings to factory defaults and all the settings will lost forever. The unit will restart at the same time.
- You can go to next parameter setting by pressing  , or pressing  key to quit and go back to measurement mode.

6.4 Setting Parameter




1. Long press  to enter into Setting parameter when in measurement mode. Press   to check through all the parameters.
2. Press  to quit and go back to measurement mode.






6.5 Calibration

Note: Calibration must be done in the following three occasions:

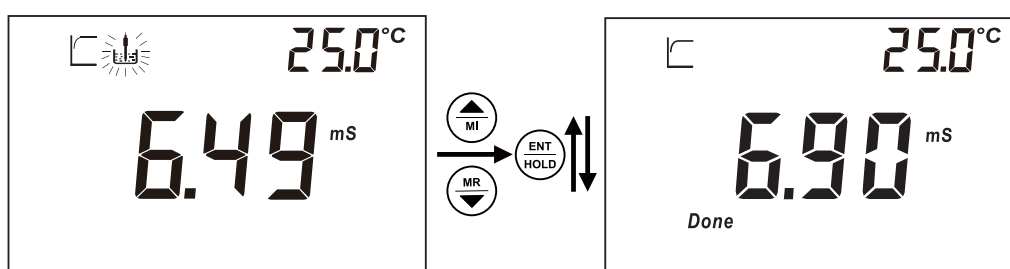1. Before using new sensors or new meter.
2. When the sensor been used for a long time and cause a wide measuring result.
3. When reference temperature or temperature factor need to be changed.

Calibration step



1. In measurement mode, press  to enter calibration mode.
2. Dip the electrode into standard solution.
3. After you get a stable current value or the  is showed, press   to adjust the value figures to match with the standard solution. Adjustable range is +30% of the stable current value. Press  to finish the calibration process.
4. The value of cell constant K indicates the status of the electrode.

6.6 Appendix

6.6.1 Conductivity sensor measuring range:

Cell Constant	Range	Resolution
k=0.01 cm	0.000~4.000 $\mu\text{S}/\text{cm}$	0.001 $\mu\text{S}/\text{cm}$
	0.00~40.00 $\mu\text{S}/\text{cm}$	0.01 $\mu\text{S}/\text{cm}$
	0.0~400.0 $\mu\text{S}/\text{cm}$	0.1 $\mu\text{S}/\text{cm}$
k=0.1 cm	0.00~40.00 $\mu\text{S}/\text{cm}$	0.01 $\mu\text{S}/\text{cm}$
	0.0~400.0 $\mu\text{S}/\text{cm}$	0.1 $\mu\text{S}/\text{cm}$
	0~4000 $\mu\text{S}/\text{cm}$	1 $\mu\text{S}/\text{cm}$
k=1 cm	0.0~400.0 $\mu\text{S}/\text{cm}$	0.1 $\mu\text{S}/\text{cm}$
	0~4000 $\mu\text{S}/\text{cm}$	1 $\mu\text{S}/\text{cm}$
	0~40.00 mS/cm	0.01 mS/cm
k=10 cm	0~4000 $\mu\text{S}/\text{cm}$	1 $\mu\text{S}/\text{cm}$
	0~40.00 mS/cm	0.01 mS/cm
	0~400.0 mS/cm	0.1 mS/cm

6.6.2 Factory Defaults

Instrument Constant	Measurement Range	Initial value
Cell constant type	0.01, 0.1, 1, 10	1.0
Temperature Base	15.0~35.0 °C	25.0 °C
Conductivity variance ratio per temperature	0.00~10.00%	2.00%
TDS factor	0.40~1.00	0.50
Salinity factor	0.48~0.65	0.65
Temperature bias	+10 °C of the indicated temperature	0.0 °C
Auto Lock function	OFF or ON	OFF
Auto Power off function	ON or OFF	ON

Dissolved Oxygen

DO200 Portable Meter

DO500 Benchtop Meter

2014 New Model! Designed for all-weather precise, easy operation comes with high luminant backlight. Ideal for testing water quality, wastewater, aquaculture and broadly used in fermentation, offers high accuracy and best value measurement experience.

- All-weather precise, Comfortable holding, Easy carrying , Simple Operation
- 65*40mm, large LCD with backlight for easy meter information reading
- IP67 rated, dustproof and waterproof, Floats on water - **[DO200]**
- Unit display: mg/L or %
- One key to check through all the settings, including: Zero offset, slope and all the settings
- Galvanic cell mode, without polarization, easy measuring
- Auto lock function
- 50 data sets memory stores and recalls
- Auto Power off saves battery after 10-mins non-use
- 2*1.5V 7AAA battery, long battery life - **[DO200]**
- CP337 Operating Carrying Pouch - **[DO200]**
- Economical, easy to operate, cost saving
- Detachable Electrode Stand organizes multiple electrodes neatly, easy installing at either left or right side and holds them firmly in place - **[DO500]**

7.1 Technical Specifications

		DO200	DO500
Testing items	Range	0.0% ~ 400.0%	
	Resolution	0.1%	
	Accuracy	±0.2% FS	
Oxygen concentration	Range	0.00 mg/L ~ 40.00 mg/L	
	Resolution	0.01 mg/L	
	Accuracy	±0.2% FS	
Temperature	Range	when compensating: 0.0 °C ~ 50.0 °C when testing: 0.0 °C ~ 50.0 °C	
	Resolution	0.1 °C	
	Accuracy	±0.2 °C	
Settings	Range	600 mbar ~ 1400 mbar	
Atmospheric Pressure	Resolution	1 mbar	
	Default	1013 mbar	
Salinity	Range	0.0 g/L ~ 40.0 g/L	
	Resolution	0.1 g/L	
	Default	0.0 g/L	
Others	Power supply	2*1.5V 7AAA batteries	Transformer 100 ~ 240V
	Working time	>500 hrs	—
Dimensions	Environment	-5 °C ~ 60 °C , Relative humidity<90%	
	Memory	50 sets	
Weight		94*190*35mm (W*L*H)	140*210*35mm (W*L*H)
		250g	650g

7.2 Browse the parameter


Long press  for 3 seconds in measurement mode.

Then press  or  to browse the parameters.

1. Working voltage
2. Slope
3. Zero offset
4. Manual Temperature compensation
5. Atmospheric Pressure
6. Salinity
7. Hold Auto Lock
8. Auto, power off
9. Battery power - **[DO200]**

7.3 Set Up

7.3.1 P01 Atmospheric Pressure

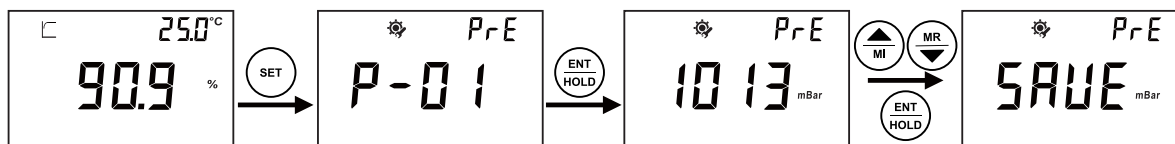




- In measurement mode, press to switch between Saturation Percentage and Oxygen concentration mg/L. Press to enter into parameter setting mode. Press to enter into electrode calibration mode.
- In P-01, you can adjust Atmospheric Pressure value for compensation to get the better measurement result.
- Please refer to above description steps to set up P-01.
- Range: 600mbar-1400mbar.
- You can go to next parameter setting by pressing , or press key to quit and go back to measurement mode.

Note:

1. During setting, you can press key to quit setting whenever necessary.
2. During setting, you can adjust anything twinkling by pressing . If it is the data twinkling, you can speedy adjust the data by pressing .




7.3.2 P02 Salinity Compensation




- In P-02, you can set up Salinity concentration value for compensation to get the better measurement result.
- Please refer to above description steps to set up P-02.
- Range: 0.0-40.0mg/L.
- You can go to next parameter setting by pressing   , or press  key to quit and go back to measurement mode.






7.3.3 P03 Temperature Offset

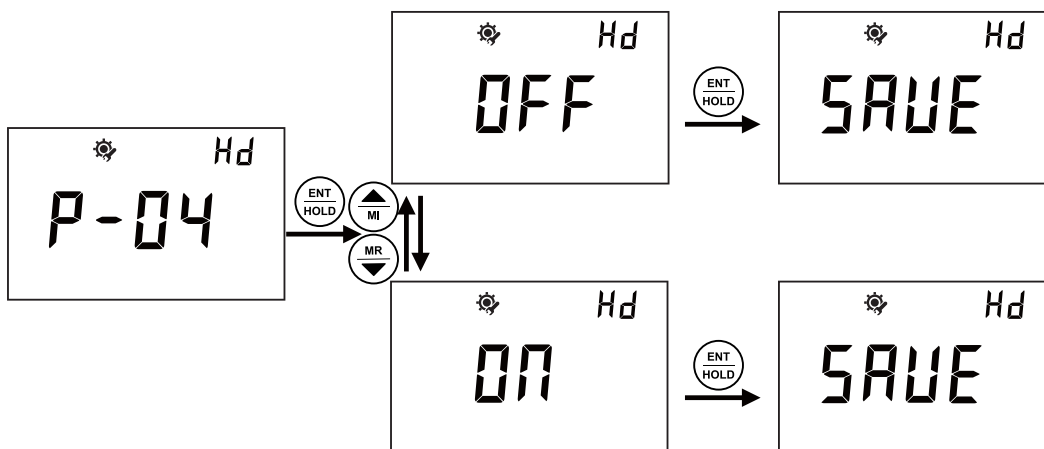


- In P-03, you can set up temperature offset value for better measurement result.
- Please refer to above description steps to set up P-03.
- Range: +10°C.
- You can go to next parameter setting by pressing   , or press  key to quit and go back to measurement mode.




7.3.4 P04 Auto Lock Function Setting




- In P-04, you can set the Auto Lock Function.
- Please refer to above description steps to set up P-04.
- If the Auto Lock Function is ON, the meter can lock the measured value after the reading had stabilized and display . Press  to unlock.
- You can go to next parameter setting by pressing   , or press  key to quit and go back to measurement mode.

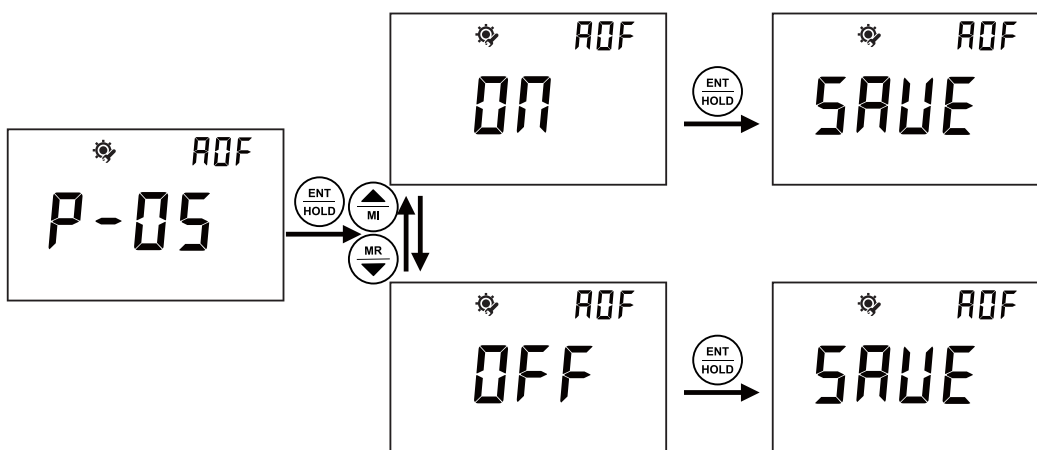7.3.5 P05 Auto Power Off Setting




- In P-05, you can set the Auto Power Off Function.
- Please refer to above description steps to set up P-05.
- If the Auto Power Off is ON, the meter will power off if no operation in 10mins.
- You can go to next parameter setting by pressing   , or press  key to quit and go back to measurement mode.

7.3.6 P06 Clearing Memory Function

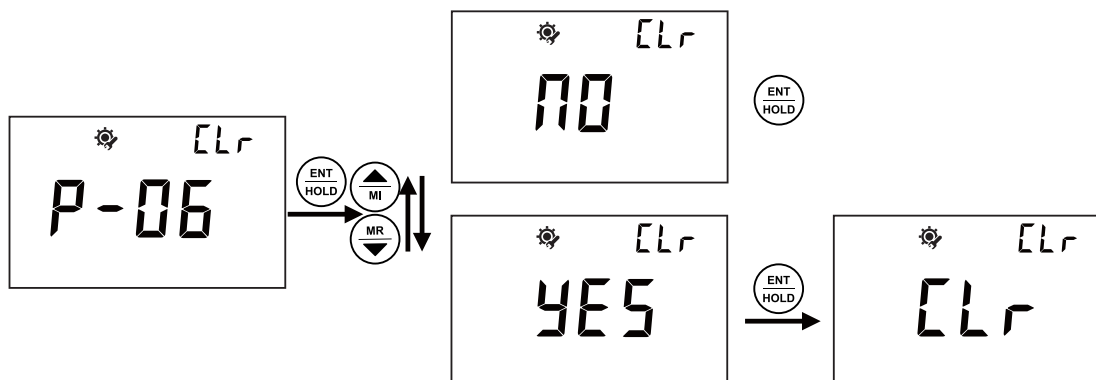


- In P-06, you can set the Clearing Memory Function.
- Please refer to above description steps to set up P-06.
- If select YES, the meter will clear all the saved data.
- You can go to next parameter setting by pressing , or press key to quit and go back to measurement mode.

7.3.7 P07 Reverting to Factory Default Setting

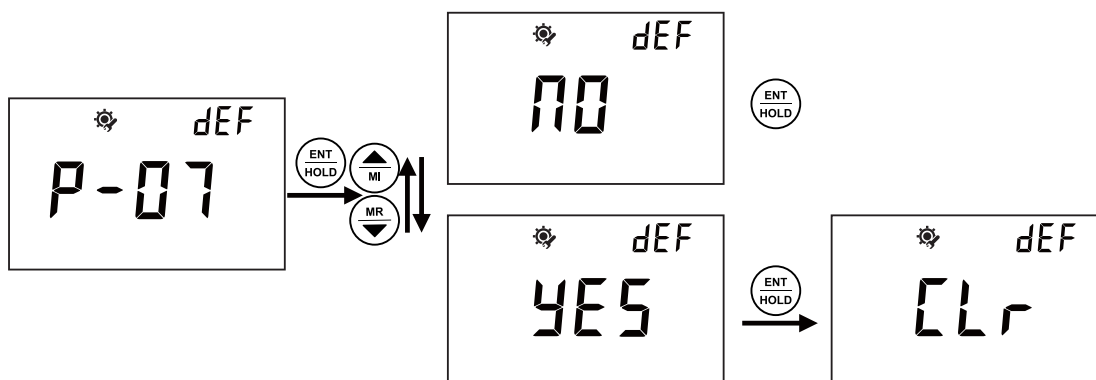

- In P-07, you can select NO or YES to not reverting or reverting to factory settings by pressing.
- Please refer to above description steps to set up P-07.
- If select YES, the meter will revert all the settings to factory defaults and all the settings will lost forever. The unit will restart at the same time.
- You can go to next parameter setting by pressing , or pressing key to quit and go back to measurement mode.

7.4 Calibration

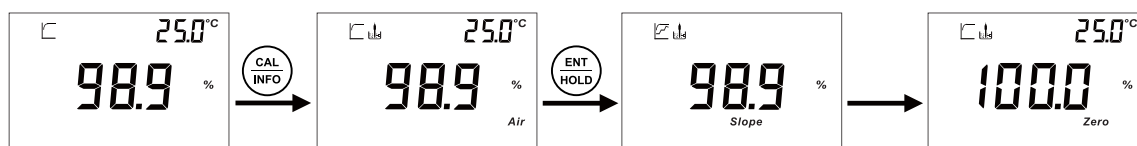

1. In measurement mode, press **CAL INFO** to enter calibration process.
2. First point calibration-100% air calibration: put the sensor in air, after you get a stable value or the stable symbol shows up, press **ENT HOLD** key to confirm 100.0% saturated dissolved oxygen point.
3. After pressing **ENT HOLD**, the electrode Slope will be showed on the screen. Then automatically initiate second point calibration -0% point calibration.

Note: Factory defaults first point 100% air calibration. If you don't need to calibrate 0% point, press **MODE ESC** key to go back to measurement mode.

After air calibration, the meter will automatically initiate second point calibration:



1. Dip the sensor in standard solution, when the value is stable, press **ENT HOLD** key to confirm 0% point. Then system go back to measurement mode.
2. If you need to adjust the Atmospheric Pressure Compensation and the Salinity Compensation according to local situation. You can set in P-01 and P-02.





Illustration on the calibration point symbols:

1. The "Air" is shown at the right bottom of the screen during and after air calibration.
2. The "Zero" is shown at the right bottom of the screen during and after air calibration.
3. Both "Air" and "Zero" is shown after 2 points calibration.

Note:

You can press **MODE ESC** back to the measurement mode during calibration and all the calibration point is saved. When the measured signal is out-ranged, the sign "Err" will be showed, possibly the aging electrode or the polluted standard solutions cause this.




7.5 Measurement Mode

1. You can first set the meter per your request. Otherwise, the setting will be factory defaults.
2. Before each measuring, first lay aside the electrode for around 15mins to fully activate the sensor after the electrode is connected with the meter.
3. After power on, press  to select the operation mode (% or mg/L).
4. Immerse the electrode in the solution and stir gently. The measuring mark will twinkle. You can read the value once the data is stabilized and shows the stable symbol .
5. If the Auto Lock Function is ON, the meter can lock the measured value after the reading had stabilized and display . Press  to unlock.
6. If the changed temperature is needed, you can refer to P-03 Temperature Offset.

7.6 Storage of Data


7.6.1 Storage of measuring data







1. If you need to save the measuring data once it is stabilized, press  , the Lot no. will be shown on the upper right of the screen. Press " " to save the data. You can also press  to not save. The meter will automatically go back to measurement mode after data saved. The unit at most can save 50 sets data. The screen will show "OVER" if more than 50 sets reminds you the memory card is full.

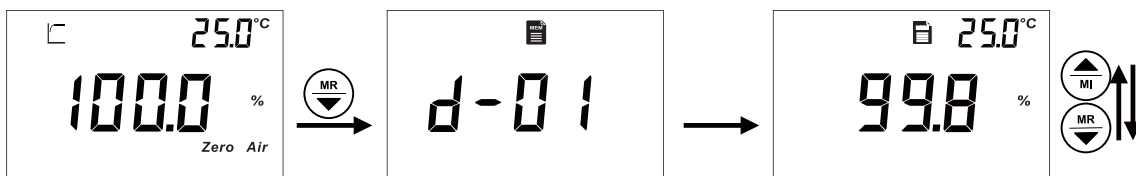2. Enter into setting mode P-06 if you need to delete the saved data forever. You can refer to Clearing memory function.


7.6.2 Browse on data storage



1. Press  to enter into the browse on the data storage , the Lot no. will be shown and data twinkling. Press   to check all the saved data. Press  to go back to measuring mode. If showed "NONE" means no data is saved yet.

2. If you quit and re-entering during checking the data, the meter will freeze in the last set of data.

7.6.3 Setting parameter

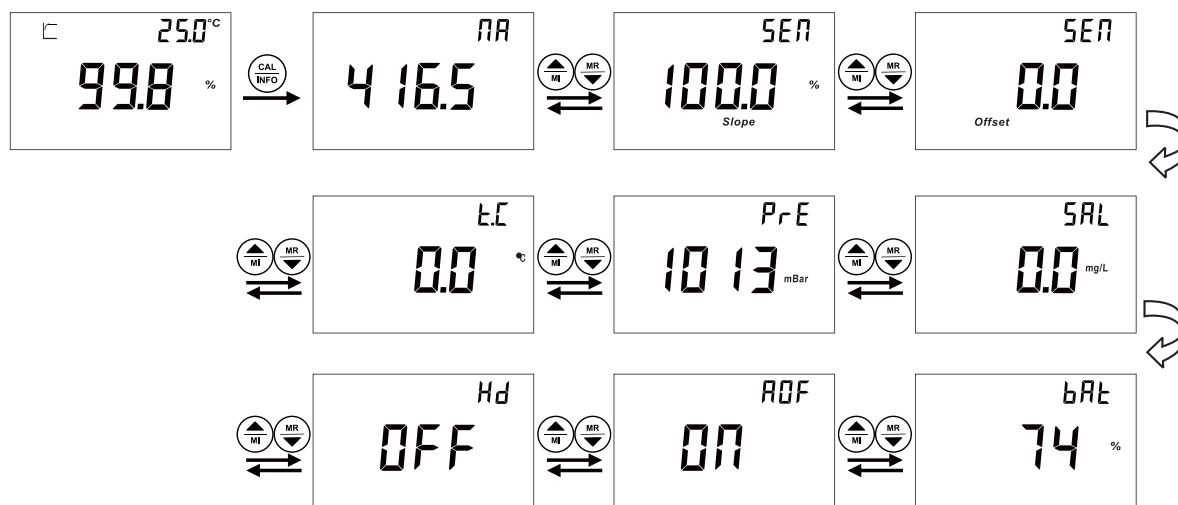

1. Long press to enter into Setting parameter when in measurement mode. Press to check through all the parameters.
2. Press to quit and go back to measurement mode.

7.7 APPENDIX

7.7.1 The corresponding Average Atmospheric Pressure with the Altitude

Altitude h/m	Average Atmospheric Pressure ph/kPa	Altitude h/m	Average Atmospheric Pressure ph/kPa
0	101.3	2800	71.4
100	100.1	2900	70.5
200	98.8	3000	69.6
300	97.6	3100	68.7
400	96.4	3200	67.9
500	95.2	3300	67.0
600	94.0	3400	66.2
700	92.8	3500	65.4
800	91.7	3600	64.6
900	90.5	3700	64.6
1000	89.4	3800	63.0
1100	88.3	3900	62.2
1200	87.2	4000	61.4
1300	86.1	4100	60.7
1400	85.0	4200	59.9
1500	84.0	4300	59.2
1600	82.9	4400	58.4
1700	81.9	4500	57.7
1800	80.9	4600	57.0
1900	79.9	4700	56.3
2000	78.9	4800	55.6
2100	77.9	4900	54.9
2200	76.9	5000	54.2
2300	76.0	5100	53.5
2400	75.0	5200	52.9
2500	74.1	5300	52.2
2600	73.2	5400	51.6
2700	72.3	5500	50.9

7.7.2 Water saturated dissolved oxygen in different temperature

(Atmospheric Pressure 101325Pa 100% RH Salinity 0‰)

Temperature°C	Dissolved oxygen(mg/L)	Temperature°C	Dissolved oxygen(mg/L)
0	14.64	20	9.08
1	14.22	21	8.90
2	13.82	22	8.73
3	13.44	23	8.57
4	13.09	24	8.41
5	12.74	25	8.25
6	12.42	26	8.11
7	12.11	27	7.96
8	11.81	28	7.82
9	11.53	29	7.69
10	11.26	30	7.56
11	11.01	31	7.43
12	10.77	32	7.30
13	10.53	33	7.18
14	10.30	34	7.07
15	10.08	35	6.95
16	9.86	36	6.84
17	9.66	37	6.73
18	9.46	38	6.63
19	9.27	39	6.53

7.7.3 Factory defaults

Item	Corresponding interface	Factory Defaults
Atmospheric Pressure	P - 01	1013 mBar
Salinity	P - 02	0.0 g/L
Auto Lock	P - 04	OFF
Auto power off	P - 05	ON
Backlight		OFF
Unit of the displayed value		%

8 GENERAL INFORMATION

8.1 Warranty

CLEAN Instruments warrants this product to be free from significant deviations in material and workmanship for a period of one year from the date of purchase. If repair is necessary and has not been the result of abuse or misuse within the warranty period, please return to CLEAN Instruments and amendment will be made without any charge. CLEAN Instruments Customer Service Center will determine if product problem is due to deviations or customer abuse. Out of warranty products will be repaired on a charge basis.

8.2 Return Of Malfunction Instruments

Authorization must be obtained from CLEAN Instruments Customer Service Center to issue a RIR number before returning items for any reason. When applying for authorization, please include date requiring the reason of return. Instruments must be carefully packed to prevent damage in shipment and insured against possible damage or loss. CLEAN Instruments will not be responsible for any damage resulting from careless or insufficient packing.

Warning: Damage as a result of inadequate packaging is the User / distributor's responsibility. Please follow the guidelines below before transporting.

8.3 Guidelines Or Returning Unit For Repair

Use the original packaging material if possible, when transporting back the unit for repair. Otherwise wrap it with bubble pack and use a corrugated box for better protection. Include a brief description of any faults suspected for the convenience of Customer Service Center, if possible. If there are any questions, feel free to contact our Customer Service Center or distributors.

*** For more new products, please feel free to contact our local distributors or log in our website.**



PH200/CON200/DO200 Portable Meter
PH500/CON500/DO500 Benchtop Meter

www.cleaninst.com

E-mail: CS@cleaninst.com

CLEAN INSTRUMENTS

2006, No.511 Tianmu W. Rd. Shanghai 200070, China
2F No.140, Zhongxiao St., Zhonghe Dist., New Taipei City ,Taiwan



OStriders